

QIAEX II Handbook

For
DNA Extraction from Agarose Gels
DNA Extraction from Polyacrylamide Gels
Desalting and Concentrating DNA from Solutions

February 1999



Trademarks

Patented or patent-pending technology and /or registered or registration-pending trademarks of QIAGEN:
QIAGEN[®], QIAEX[®], QIAquick[™].

The PCR process is covered by U.S. patents 4,683,195 and 4,683,202 and foreign equivalents owned by Hoffman-La Roche AG.

Whatman is a registered trademark of Whatman Ltd.

© 1999 QIAGEN, all rights reserved.

Germany	QIAGEN GmbH	Max-Volmer-Straße 4 • 40724 Hilden Orders 02103-892-230 • Fax 02103-892-233 • Technical 02103-892-240
USA	QIAGEN Inc.	28159 Avenue Stanford • Valencia • CA 91355 Orders 800-426-8157 • Fax 800-718-2056 • Technical 800-DNA-PREP (800-362-7737)
Australia	QIAGEN Pty Ltd ACN 072 382 944	PO Box 25 • Clifton Hill • Victoria 3068 Orders 03-9489-3666 • Fax 03-9489-3888 • Technical 03-9489-3666
Canada	QIAGEN Inc.	2900 Argentia Road • Unit 23 • Mississauga • Ontario • L5N 7X9 Orders 800-572-9613 • Fax 800-713-5951 • Technical 800-DNA-PREP (800-362-7737)
France	QIAGEN S. A.	3 avenue du Canada • LP 809 • 91974 Courtaboeuf Cedex Orders 01-60-920-920 • Fax 01-60-920-925 • Technical 01-60-920-930
Japan	QIAGEN K.K.	Hakusan Takayanagi Bldg. 8F • 7-6, Hakusan 1 Chome • Bunkyo-ku, Tokyo 113-0001 Telephone 03-5805-7261 • Fax 03-5805-7263 • Technical 03-5805-7261
Switzerland	QIAGEN AG	Auf dem Wolf 39 • 4052 Basel Orders 061-319-30-30 • Fax 061-319-30-33 • Technical 061-319-30-31
UK	QIAGEN Ltd.	Boundary Court • Gatwick Road • Crawley • West Sussex, RH10 2AX Orders 01293-422-911 • Fax 01293-422-922 • Technical 01293-422-999

<http://www.qiagen.com>

Contents

Kit Contents	4
Storage Conditions	4
Product Use Limitations	4
Product Warranty and Satisfaction Guarantee	5
Quality Control	5
Technical Assistance	5
Introduction	6
About this Handbook	6
QIAEX II Specifications	8
The QIAEX II Principle	8
Solubilization of agarose without NaI	8
Efficient DNA adsorption to QIAEX II — salt and pH dependence	9
pH indicator in solubilization & binding Buffer QX1	10
Washing	10
Elution in low-salt solutions	11
Applications using QIAEX II purified fragments	11
QIAEX II Protocols	12
■ QIAEX II Agarose Gel Extraction Protocol	12
■ QIAEX II Polyacrylamide Gel Extraction Protocol	14
■ QIAEX II Protocol for Desalting and Concentrating DNA Solutions	16
Troubleshooting Guide	18
Ordering Information	21
International Sales and Distributors	23

Kit Contents

QIAEX II Gel Extraction Kit	QIAEX II (150)	QIAEX II (500)
Catalog No.	20021	20051
No. of preparations* (10 µl QIAEX II per prep)	150	500
QIAEX II Suspension	3 x 0.5 ml	5 x 1 ml
Buffer QX1 [†] (with pH indicator)	2 x 100 ml	6 x 100 ml
Buffer PE (concentrate)	2 x 20 ml	6 x 20 ml
Handbook	1	1

* Routine purifications from gel slices >100 mg which contain fragments <100 bp or where the gels are >2% agarose require additional Buffer QX1 to perform the full number of extractions (see ordering information, page 21).

[†] Buffer QX1 contains chaotropic salt which is an irritant. Take appropriate laboratory safety precautions and wear gloves when handling.

Storage Conditions

QIAEX[®] II Gel Extraction Kits should be stored dry at room temperature (15–25°C). When stored under these conditions and handled appropriately, the kits can be stored for at least 12 months without showing any reduction in performance, capacity, or quality of separation.

Product Use Limitations

QIAEX II Gel Extraction Kits are developed, designed, and sold for research purposes only. They are not to be used for human diagnostic or drug purposes or to be administered to humans unless expressly cleared for that purpose by the Food and Drug Administration in the USA or the appropriate regulatory authorities in the country of use. All due care and attention should be exercised in the handling of many of the materials described in this text.

Product Warranty and Satisfaction Guarantee

QIAGEN guarantees the performance of all products in the manner described in our product literature. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, QIAGEN will replace it free of charge or refund the purchase price. We reserve the right to change, alter, or modify any product to enhance its performance and design. If a QIAGEN product does not meet your expectations, simply call your local Technical Service Department. We will credit your account or exchange the product — as you wish.

A copy of QIAGEN terms and conditions can be obtained on request, and is also provided on the back of our invoices. If you have questions about product specifications or performance, please call QIAGEN Technical Services or your local distributor.

Quality Control

As part of the stringent QIAGEN quality assurance program, the performance of QIAEX II Kits is monitored routinely and on a lot-to-lot basis. QIAEX II Kits are tested by isolation of DNA fragments of various sizes from either aqueous solution or agarose gel. The quality of the isolated DNA is checked by several assays commonly used for nucleic acids. The DNA binding capacity of QIAEX II resin is tested by determining the recovery from a specific amount of loaded DNA.

Technical Assistance

At QIAGEN we pride ourselves on the quality and availability of our technical support. Our Technical Service Departments are staffed by experienced scientists with extensive practical and theoretical expertise in molecular biology and the use of QIAGEN products. They are always available to discuss any general or specific questions you may have. If you have any questions or experience any problems regarding the QIAEX II Gel Extraction Kit, or QIAGEN products in general, please do not hesitate to contact us.

QIAGEN customers are also a major source of information regarding advanced or specialized uses of our products. This information is helpful to other scientists as well as to the researchers at QIAGEN. We therefore also encourage you to contact us if you have any suggestions about product performance or new applications and techniques.

For technical assistance and more information please call one of the QIAGEN Technical Service Departments or local distributors listed on the last page.

Introduction

QIAEX II Gel Extraction Kit is designed to extract and purify DNA from any agarose gel in either TAE (Tris-acetate/EDTA) or TBE (Tris-borate/EDTA) buffer without phenol extraction or ethanol precipitation. The QIAEX II silica particles have been optimized to enhance recovery of very small and very large DNA fragments. DNA molecules of 40 bp to 50 kb are adsorbed to QIAEX II particles in the presence of high salt. All non-nucleic acid impurities such as agarose, proteins, salts, and ethidium bromide are removed during washing steps. The pure DNA is efficiently eluted in just 20 μ l of Tris buffer or water and is suitable for most subsequent applications, e.g., restriction digestion, labeling, ligation, PCR, sequencing, and microinjection (see Applications using QIAEX II Purified Fragments, page 11).

About this Handbook

This handbook contains technical information about QIAEX II to help you make the most of the system. The handbook explains the principle behind QIAEX II and outlines the major steps of the procedure. Detailed protocols for agarose gel extraction, polyacrylamide gel extraction, and for desalting and concentrating DNA solutions are provided. A troubleshooting guide is also included to help analyze problems which may occur.

QIAEX II adds a number of new advantages such as simplified handling and a more robust, streamlined procedure, to all the familiar benefits of its predecessor — QIAEX. QIAEX II silica particles are even easier to pellet and resuspend, and are also less susceptible to over-drying before elution. The QIAEX II Gel Extraction Kit also requires fewer wash steps and buffers. The solubilization and binding Buffer QX1 now contains a pH indicator, allowing easy determination of optimal pH for DNA-binding. Turn to page 10 for detailed information.

For long-time QIAEX users, please become familiar with the appropriate new protocol in this handbook and discard any old QIAEX handbooks in the laboratory.

IMPORTANT: Always use the protocol or handbook provided with the product.

QIAEX II Gel Extraction Procedure

Solubilized gel slice
+ QIAEX II particles



Bind



Collect



Wash



Elute



Pure DNA Fragment

QIAEX II Specifications

Maximum binding capacity:	5 µg DNA per 10 µl QIAEX II suspension
Recovery:	60–95% of DNA fragments 40 bp – 50 kb
Minimum elution volume:	20 µl

The QIAEX II Principle

With the QIAEX II Gel Extraction Kit, extraction and purification of DNA fragments are based on solubilization of agarose and selective, quantitative adsorption of nucleic acids to the QIAEX II silica-gel particles in the presence of high salt. Elution of the DNA is accomplished with a low-salt solution such as Tris buffer or water.

Solubilization of agarose without NaI

The optimized Buffer QX1 in QIAEX II Kits efficiently solubilizes agarose, and does not contain NaI. Residual NaI may be difficult to remove from DNA samples, and reduces the efficiency of subsequent enzymatic reactions such as blunt-end ligation.

The standard QIAEX II Gel Extraction protocol is used to extract DNA from 0.3–2% standard or low-melt agarose gels in TAE or TBE buffer. The concentration of agarose in the gel determines the size range of DNA molecules which can be resolved (Table 1).

Table 1. Separation range in TAE gels containing different concentrations of agarose (1)

Agarose	Size of DNA fragments separated
0.3%	5 – 60 kb
0.5%	1 – 30 kb
0.7%	0.8 – 12 kb
1.0%	0.5 – 10 kb
1.2%	0.4 – 7 kb
1.5%	0.2 – 3 kb
2.0%	0.05 – 2 kb

A typical agarose gel slice is solubilized by adding 3 volumes of Buffer QX1 to 1 volume of gel (e.g., 300 μ l of Buffer QX1 is added to 100 mg gel slice) and incubating at 50°C for 10 minutes. The high concentration of a chaotropic salt in Buffer QX1 disrupts hydrogen bonding between sugars in the agarose polymer, allowing solubilization of the gel slice. In addition, the high salt concentration dissociates DNA binding proteins from the DNA fragments.

With Buffer QX1, DNA can be extracted from TBE gels without an extra TBE modifier. TBE gels are not as readily solubilized as TAE gels due to the formation of complexes between borate and the cis-diol groups of sugar monomers and polymers. Chaotropic salts cannot solubilize these gels without additional sugar monomers such as sorbitol or mannitol to chelate the free borate. Ready-to-use Buffer QX1 has been optimized to effectively dissolve either TAE or TBE gels.

A \leq 2% agarose gel slice is normally solubilized within 2–3 minutes with Buffer QX1 at 50°C. The incubation with QIAEX II is extended to 10 minutes to complete the adsorption of DNA to the QIAEX II particles.

Efficient DNA adsorption to QIAEX II — salt and pH dependence

Adsorption of DNA to glass, silica gel, or diatomaceous earth in high salt is a well known phenomenon (2). Incubation of a DNA solution in a highly electrolytic environment with large anions (e.g., from chaotropic salts) causes a modification in the structure of water (3), forcing the DNA to adsorb to the silica particles. Adsorption of fragments smaller than 100 bp is enhanced by increasing the salt concentration, while fragments larger than 4 kb are adsorbed at lower salt concentrations.

Adsorption of DNA to silica also depends on pH. Adsorption efficiency is typically 95% if the pH is \leq 7.5, and is drastically reduced at higher pH (Figure 1). If the pH of the binding mixture is $>$ 7.5, the optimal pH for DNA binding can be achieved by adding a small volume of 3M sodium acetate, pH 5.0.

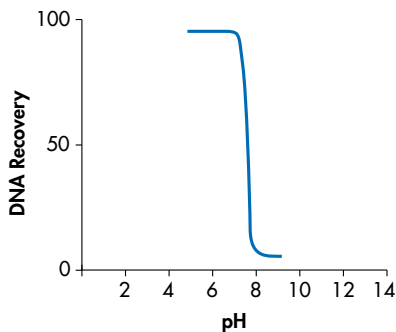


Figure 1. pH dependence of DNA adsorption to silica. 1 μ g of a 2.9-kb fragment was adsorbed at different pH values and eluted with 10 mM Tris-Cl, pH 8.5. The graph shows percentage DNA recovery, reflecting the relative adsorption efficiency, versus pH of adsorption.

pH indicator in solubilization & binding Buffer QX1

Buffer QX1 in QIAEX II Gel Extraction Kits is used for solubilization of agarose gel slices and binding of DNA to the QIAEX II silica particles. Buffer QX1 now contains a pH indicator, allowing easy determination of the optimal pH for DNA binding. DNA adsorption requires a pH ≤ 7.5 , and the pH indicator appears yellow in this range. If the pH is >7.5 , which can occur if the agarose gel electrophoresis buffer is frequently used or incorrectly prepared, the binding mixture turns orange or violet (Figure 2). This means that the pH of the sample exceeds the buffering capacity of Buffer QX1 and DNA adsorption will be inefficient. In this case, the pH of the binding mixture can easily be corrected by addition of a small volume of 3M sodium acetate, pH 5.0, before proceeding with the protocol.

As an additional benefit, the color of the binding mixture allows easy visualization of any unsolubilized agarose, ensuring complete solubilization and maximum yields. The indicator dye does not interfere with DNA binding and is completely removed during the cleanup procedure.

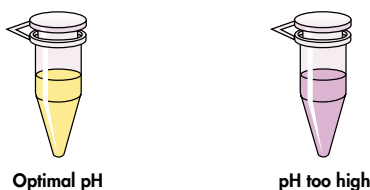


Figure 2. Indicator dye in solubilization & binding Buffer QX1 identifies optimal pH for DNA binding.

Washing

DNA molecules bind to the QIAEX II particles during the adsorption step, and all non-nucleic acid impurities such as agarose, proteins, ethidium bromide, and salts remain in the supernatant. A high salt wash with Buffer QX1 removes residual agarose, and two washes with ethanol-containing Buffer PE efficiently remove salt contaminants. All traces of supernatant must be carefully removed at each step to eliminate impurities and reduce buffer carryover.

Washed QIAEX II particles carrying adsorbed DNA are pelleted and air-dried at room temperature for 10–15 minutes. Drying the pellet is necessary to remove all traces of residual ethanol which may interfere with subsequent enzymatic reactions.

Elution in low-salt solutions

Elution efficiency depends on pH and salt concentration. Contrary to adsorption, elution is best under basic conditions and low salt concentrations when using QIAEX II. DNA is typically eluted with 20 μ l of 10 mM Tris-Cl, pH 8.5, or with water. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water to elute, make sure that the pH value is within this range. In addition, DNA must be stored at -20°C when eluted with water, since DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE (10 mM Tris-Cl, 1 mM EDTA, pH 8.0), but the EDTA may inhibit subsequent enzymatic reactions.

Elution efficiency also depends on temperature. DNA fragments smaller than 4 kb are recovered with 70–95% efficiency when eluted at room temperature with a 5-min incubation. For efficient elution of DNA fragments larger than 4 kb, the incubation temperature should be increased to 50°C .

After elution, QIAEX II is pelleted by centrifugation. The supernatant containing pure DNA is easily removed from the QIAEX II pellet, ready for use in subsequent applications.

Applications using QIAEX II purified fragments

DNA purified with QIAEX II is suitable for various applications, for example, restriction digestion, labeling, ligation, transformation, PCR, sequencing, *in vitro* transcription, and microinjection. Some applications, such as sequencing, blunt-end ligation, *in vitro* transcription, or microinjection, are highly sensitive to resin carryover and residual salts, and therefore all traces of supernatant must be carefully removed at each wash step. After elution, the supernatant containing the pure DNA should be carefully removed from the QIAEX II pellet. If necessary, the supernatant may be centrifuged a second time to ensure complete removal of all QIAEX II particles.

For sensitive downstream applications such as fluorescent sequencing, QIAquick™ Gel Extraction Kits (see ordering information, page 21) are recommended. QIAquick Gel Extraction Kits combine a convenient microspin format with a specially designed silica membrane. The quality of the purified DNA is less dependent on the handling procedures.

1. Sambrook, J., Fritsch, E.F., and Maniatis, T. (1989) *Molecular Cloning: A Laboratory Manual*, Cold Spring Harbor, NY: Cold Spring Harbor Laboratory Press.
2. Vogelstein, B., and Gillespie, D. (1979) Preparative and analytical purification of DNA from agarose. *Proc. Natl. Acad. Sci. USA* **76**, 615–619.
3. Hamaguchi, K., and Geiduschek, E.P. (1962) The effect of electrolytes on the stability of deoxyribonucleate helix. *J. Am. Chem. Soc.* **84**, 1329–1337.

QIAEX II Agarose Gel Extraction Protocol

This protocol is designed for the extraction of 40-bp to 50-kb DNA fragments from 0.3–2% standard or low-melt agarose gels in TAE or TBE buffers.

- Notes:**
- The yellow color of Buffer QX1 indicates a pH ≤ 7.5 .
 - Add ethanol (96–100%) to Buffer PE before use (see bottle label for volume).
 - A heating block or water bath at 50°C is required.
 - 3M sodium acetate, pH 5.0, may be necessary.
 - All centrifugation steps are at maximum speed ($\geq 10,000 \times g$, $\sim 13,000$ rpm) in a conventional, table-top microcentrifuge.
 - For DNA fragments larger than 10 kb, mix by gently flicking the tube to avoid shearing the DNA. Do not vortex the tube.

1. Excise the DNA band from the agarose gel with a clean, sharp scalpel.

Minimize the size of the gel slice by removing excess agarose. Use a 1.5-ml microfuge tube for processing up to 250 mg agarose.

2. Weigh the gel slice in a colorless tube. Add 3 volumes of Buffer QX1 to 1 volume of gel for DNA fragments 100 bp – 4 kb; otherwise, follow the table below.

For example, add 300 μ l of Buffer QX1 to each 100 mg of gel.

DNA fragments <100 bp	Add 6 volumes of Buffer QX1
DNA fragments >4 kb	Add 3 volumes of Buffer QX1 plus 2 volumes of H ₂ O
>2% or Metaphor agarose gels	Add 6 volumes of Buffer QX1

3. Resuspend QIAEX II by vortexing for 30 sec. Add QIAEX II to the sample according to the table below and mix.

≤ 2 μ g DNA	Add 10 μ l of QIAEX II
2–10 μ g DNA	Add 30 μ l of QIAEX II
Each additional 10 μ g DNA	Add additional 30 μ l of QIAEX II

4. Incubate at 50°C for 10 min to solubilize the agarose and bind the DNA. Mix by vortexing* every 2 min to keep QIAEX II in suspension. Check that the color of the mixture is yellow.

If the color of the mixture is orange or purple, add 10 μ l 3M sodium acetate, pH 5.0, and mix. The color should turn to yellow. The incubation should then be continued for an additional 5 min at least.

The adsorption of DNA to QIAEX II particles is only efficient at pH ≤ 7.5 . Buffer QX1 now contains a pH indicator which is yellow at pH ≤ 7.5 , and orange or violet at higher pH, allowing easy determination of the optimal pH for DNA binding.

5. Centrifuge the sample for 30 sec and carefully remove supernatant with a pipet.

6. Wash the pellet with 500 μ l of Buffer QX1.

Resuspend the pellet by vortexing*. Centrifuge the sample for 30 sec and remove all traces of supernatant with a pipet. This wash step removes residual agarose contaminants.

7. Wash the pellet twice with 500 μ l of Buffer PE.

Resuspend the pellet by vortexing*. Centrifuge the sample for 30 sec and carefully remove all traces of supernatant with a pipet. These washing steps remove residual salt contaminants.

8. Air-dry the pellet for 10–15 min or until the pellet becomes white.

If 30 μ l of QIAEX II suspension is used, air-dry the pellet for approximately 30 min. Do not vacuum dry, as this may cause overdrying. Overdrying the QIAEX II pellet may result in decreased elution efficiency.

9. To elute DNA, add 20 μ l of 10 mM Tris-Cl, pH 8.5 or H₂O and resuspend the pellet by vortexing*. Incubate according to the table below.

DNA fragments ≤ 4 kb	Incubate at room temp. for 5 min
DNA fragments 4–10 kb	Incubate at 50°C for 5 min
DNA fragments > 10 kb	Incubate at 50°C for 10 min

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water for elution, make sure that the pH is within this range, and store DNA at -20°C as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE buffer (10 mM Tris-Cl, 1 mM EDTA, pH 8.0) but the EDTA may inhibit subsequent enzymatic reactions.

10. Centrifuge for 30 sec. Carefully pipet the supernatant into a clean tube.

The supernatant now contains the purified DNA.

11. Optional: repeat steps 9 and 10 and combine the eluates.

A second elution step will increase the yield by approximately 10–15%.

* For fragments larger than 10 kb, resuspend the pellet by inverting and flicking the tube. Vortexing can cause shearing of large DNA fragments.

QIAEX II Polyacrylamide Gel Extraction Protocol

- Notes:**
- The yellow color of Buffer QX1 indicates a pH ≤ 7.5 .
 - Add ethanol (96–100%) to Buffer PE before use (see bottle label for volume).
 - A heating block or water bath at 50°C is required.
 - 3M sodium acetate, pH 5.0, may be necessary.
 - All centrifugation steps are at maximum speed ($\geq 10,000 \times g$, ~13,000 rpm) in a conventional, table-top microcentrifuge.
 - Prepare diffusion buffer: 0.5 M ammonium acetate; 10 mM magnesium acetate; 1 mM EDTA, pH 8.0; 0.1% SDS.
 - A disposable plastic column or a syringe barrel containing either a Whatman® GF/C filter or packed, siliconized glass wool will be required for each extraction.

- 1. Excise the gel slice containing the DNA band with a clean, sharp scalpel.**
Minimize the size of the gel slice by removing excess polyacrylamide.
- 2. Weigh the gel slice. Add 1–2 volumes of diffusion buffer to 1 volume of gel (i.e., 100–200 μ l for each 100 mg of gel).**
- 3. Incubate at 50°C for 30 min.**
- 4. Microcentrifuge the sample for 1 min.**
- 5. Carefully remove the supernatant using a pipet or a drawn-out Pasteur pipet. Pass the supernatant through a disposable plastic column or a syringe containing either a Whatman GF/C filter or packed, siliconized glass wool to remove any residual polyacrylamide.**
- 6. Calculate the approximate volume of the recovered supernatant.**
- 7. For DNA fragments <100 bp, add 6 volumes of Buffer QX1 to 1 volume of sample. Otherwise, add 3 volumes Buffer QX1. Check that the color of the mixture is yellow.** If the color of the mixture is orange or purple, add 10 μ l 3M sodium acetate, pH 5.0, and mix. The color should turn to yellow.

The adsorption of DNA to QIAEX II particles is only efficient at pH ≤ 7.5 . Buffer QX1 now contains a pH indicator which is yellow at pH ≤ 7.5 , and orange or violet at higher pH, allowing easy determination of optimal pH for DNA-binding.

- 8. Resuspend QIAEX II by vortexing for 30 sec.**

9. Add 10 μ l of QIAEX II and mix. Incubate at room temperature for 10 min. Vortex every 2 min to keep QIAEX II in suspension.

10. Centrifuge the sample for 30 sec and remove supernatant.

11. Wash the pellet twice with 500 μ l of Buffer PE.

12. Air-dry the pellet for 10–15 min or until the pellet becomes white.

Do not vacuum dry, as this may cause overdrying. Overdrying the QIAEX II pellet may result in decreased elution efficiency.

13. To elute DNA, add 20 μ l of 10 mM Tris-Cl, pH 8.5, or H₂O and resuspend the pellet by vortexing. Incubate for 5 min at room temperature.

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water for elution, make sure that the pH is within this range, and store DNA at -20°C as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE buffer (10 mM Tris-Cl, 1 mM EDTA, pH 8.0) but the EDTA may inhibit subsequent enzymatic reactions.

14. Centrifuge for 30 sec. Carefully transfer the supernatant into a clean tube.

The supernatant now contains pure DNA.

15. **Optional: repeat steps 13 and 14 and combine the eluates.**

A second elution step will increase the yield by approximately 10–15%.

QIAEX II Protocol for Desalting and Concentrating DNA Solutions

QIAEX II Gel Extraction Kit can be used to purify and concentrate DNA fragments from 40 bp to 50 kb from aqueous solutions without phenol extraction or ethanol precipitation. Enzymes such as phosphatases, restriction endonucleases, polymerases, as well as dNTPs, and salts are removed. For direct purification from PCR reactions, use the QIAquick PCR Purification Kit (see ordering information, page 21) to recover PCR products and efficiently remove primers.

- Notes:**
- The yellow color of Buffer QX1 indicates a pH ≤ 7.5 .
 - Add ethanol (96–100%) to Buffer PE before use (see bottle label for volume).
 - All centrifugation steps are at maximum speed ($\geq 10,000 \times g$, $\sim 13,000$ rpm) in a conventional table-top microcentrifuge.
 - For DNA fragments larger than 10 kb, mix by gently flicking the tube to avoid shearing the DNA. Do not vortex the tube.

1. Transfer the sample to a colorless tube. Add 3 volumes of Buffer QX1 to 1 volume of sample.

For DNA fragments 100 bp – 4 kb; otherwise follow the table below.

DNA fragments <100 bp	Add 6 volumes of Buffer QX1
DNA fragments >4 kb	Add 3 volumes of Buffer QX1 plus 2 volumes of H ₂ O

2. Check that the color of the sample mixture is yellow.

If the color of the mixture is orange or purple, add 10 μ l 3M sodium acetate, pH 5.0, and mix. The color should now be yellow.

The adsorption of DNA to QIAEX II particles is only efficient at pH ≤ 7.5 . Buffer QX1 now contains a pH indicator which is yellow at pH ≤ 7.5 , and orange or violet at higher pH, allowing easy determination of the optimal pH for DNA-binding.

3. Resuspend QIAEX II by vortexing 30 sec.

4. Add 10 μ l of QIAEX II per 5 μ g of DNA and mix*. Incubate at room temperature for 10 min. Mix every 2 min to keep QIAEX II in suspension.

5. Centrifuge the sample for 30 sec and remove supernatant.

6. Wash the pellet twice with 500 μ l of Buffer PE.

7. Air-dry the pellet for 10–15 min or until the pellet becomes white.

Do not vacuum dry, as this may cause overdrying. Overdrying the QIAEX II pellet may result in decreased elution efficiency.

8. To elute DNA, add 20 μ l of 10 mM Tris-Cl, pH 8.5, or H₂O and resuspend the pellet by vortexing*. Incubate according to the table below.

DNA fragments \leq 4 kb	Incubate at room temp. for 5 min
DNA fragments 4–10 kb	Incubate at 50°C for 5 min
DNA fragments >10 kb	Incubate at 50°C for 10 min

Elution efficiency is dependent on pH. The maximum elution efficiency is achieved between pH 7.0 and 8.5. When using water for elution, make sure that the pH is within this range, and store DNA at -20°C as DNA may degrade in the absence of a buffering agent. The purified DNA can also be eluted in TE buffer (10 mM Tris-Cl, 1 mM EDTA, pH 8.0) but the EDTA may inhibit subsequent enzymatic reactions.

9. Centrifuge for 30 sec. Carefully transfer the supernatant into a clean tube.

The supernatant now contains pure DNA.

10. Optional: repeat steps 8 and 9 and combine the eluates.

A second elution step will increase the yield by approximately 10–15%.

* For fragments larger than 10 kb, resuspend the pellet by inverting and flicking the tube. Vortexing can cause shearing of large DNA fragments.

Troubleshooting Guide

This troubleshooting guide may be helpful in solving any problems which may arise. The scientists at QIAGEN Technical Services are always happy to answer any questions you may have about either the information and protocol(s) in this handbook or molecular biology applications (see page 23).

Comments and Suggestions

Poor DNA recovery

- | | | | | | | | |
|--|--|----------------------|----------------------------|------------------|----------------------------|----------------------------|---------------------------------------|
| a) Ethanol was not added to concentrated Buffer PE before use | Repeat procedure with correctly prepared Buffer PE. | | | | | | |
| b) pH of electrophoresis buffer too high (binding mixture turns orange or violet). | The electrophoresis buffer has been repeatedly used or incorrectly prepared, resulting in a sample pH that exceeds the buffering capacity of Buffer QX1 and leads to inefficient DNA binding. Add 10 μ l of 3M sodium acetate, pH 5.0, to the sample and mix. The color of the mixture should turn yellow indicating the correct pH for DNA binding. For binding mixtures with only small color changes (slight orange color), add the sodium acetate. | | | | | | |
| c) Inappropriate elution buffer | DNA will only be eluted in the presence of low salt buffer (e.g., 10 mM Tris-Cl, pH 8.5) or water. Check the pH and salt concentration of the elution buffer. | | | | | | |
| d) Incomplete solubilization of gel slice in Buffer QX1 | Mix the tube every 2 min during the solubilization step. Large gel slices may require a few extra minutes at 50°C to solubilize. DNA will remain in any undissolved agarose. | | | | | | |
| e) Low salt concentration during adsorption | Sufficient salt must be present during the adsorption step. Weigh the gel slice weight accurately and add the appropriate amount of Buffer QX1. | | | | | | |
| f) Insufficient amount of QIAEX II | Enough QIAEX II must be present at the adsorption step to bind DNA:
<table border="0" style="margin-left: 20px;"><tr><td>≤ 2 μg DNA</td><td>Add 10 μl of QIAEX II</td></tr><tr><td>2–10 μg DNA</td><td>Add 30 μl of QIAEX II</td></tr><tr><td>Each additional 10 μg</td><td>Add additional 30 μl of QIAEX II</td></tr></table> | ≤ 2 μ g DNA | Add 10 μ l of QIAEX II | 2–10 μ g DNA | Add 30 μ l of QIAEX II | Each additional 10 μ g | Add additional 30 μ l of QIAEX II |
| ≤ 2 μ g DNA | Add 10 μ l of QIAEX II | | | | | | |
| 2–10 μ g DNA | Add 30 μ l of QIAEX II | | | | | | |
| Each additional 10 μ g | Add additional 30 μ l of QIAEX II | | | | | | |

Comments and Suggestions

- g) Buffer QX1 incompletely removed before wash step with Buffer PE
If Buffer QX1 is not completely removed, the subsequent Buffer PE washing steps may not efficiently remove all traces of salt. The presence of high concentrations of salt may prevent complete elution of DNA from the QIAEX II particles. Additionally, residual salt may result in poor performance of the DNA in subsequent enzymatic reactions.
- h) Over-dried QIAEX II pellet
Do not use a vacuum dryer to dry the QIAEX II pellet. If pellet is dried too extensively, the DNA may irreversibly bind to the QIAEX II particles, resulting in decreased recovery.
- i) Insufficient volume of elution buffer
Increase volume of elution buffer in proportion to amount of extra QIAEX II added. An optional extra elution step will increase yield by approximately 10–15%.
- j) Incorrect volume of Buffer QX1 added for solubilization

<100 bp DNA	Add 6 volumes of Buffer QX1
100 bp – 4 kb DNA	Add 3 volumes of Buffer QX1
> 4 kb DNA	Add 3 volumes of Buffer QX1 plus 2 volumes of H ₂ O
> 2% or Metaphor agarose gels	Add 6 volumes of Buffer QX1

(see page 8 for a description of the underlying principle)

DNA does not perform well in subsequent reactions

- a) Incomplete removal of Buffer QX1 supernatant
Salt from Buffer QX1 may be carried over and inhibit subsequent enzymatic reactions. Wash twice with Buffer PE.
- b) Incomplete removal of Buffer PE supernatant
Residual ethanol may reduce elution efficiency. Carefully remove all Buffer PE, which contains ethanol, to enable pellet to dry sufficiently before elution.
- c) Incomplete removal of QIAEX II particles after elution
QIAEX II particles may bind proteins (e.g., enzymes) under low salt conditions, and reduce enzyme activity in subsequent reactions. If QIAEX II particles are inadvertently collected with the supernatant, centrifuge again to remove all traces of QIAEX II.

Comments and Suggestions

DNA contains residual ethanol

(samples float out of wells of agarose gels)

- a) QIAEX II pellet was under-dried
- Buffer PE was not completely removed by drying. Dry the DNA eluate briefly in a vacuum dryer to remove residual ethanol.

DNA appears to “smear” on an agarose gel.

- a) Sample well is over-loaded
- Smiling and smearing of DNA bands may occur when too much sample is loaded in one well. Use less DNA per well or use a larger well.
- b) Vigorous handling of DNA
- Vortexing and vigorous pipetting of fragments above 10 kb should be avoided throughout. QIAEX II can be easily resuspended by flicking the microfuge tube. Very large DNA fragments can be protected from shearing by keeping the QIAEX II pellet intact during the washing steps. The wash buffers should remain on the QIAEX II pellet for 5 min to allow complete diffusion.
- c) Eluate contains denatured ssDNA
- Denatured ssDNA appears as a smaller smeared band on an analytical gel. Use the eluted DNA to prepare the subsequent enzymatic reaction but omit the enzyme. To reanneal the ssDNA, incubate the reaction mixture at 95°C for 2 min, and allow the tube to cool slowly to room temperature. Add the enzyme and proceed as usual. Alternatively, the DNA can be eluted in 10 mM Tris buffer containing 10 mM NaCl. The salt and buffering agent promote the renaturation of DNA strands. However the salt concentration of the eluate must then be considered for subsequent applications.

Ordering Information

Product	Contents	Cat. No.
QIAEX II Gel Extraction Kit (150)	For 150 extractions*: 3 x 0.5 ml QIAEX II Suspension, Buffers	20021
QIAEX II Gel Extraction Kit (500)	For 500 extractions*: 5 x 1.0 ml QIAEX II Suspension, Buffers	20051
Individual Kit Components		
QIAEX II Suspension	1.5 ml QIAEX II Suspension	20902
Buffer QX1	500 ml Solubilization and Binding Buffer (with pH indicator)	20912
Buffer PE	100 ml Wash Buffer (concentrate; final volume 500 ml)	19065
Related Products		
QIAquick PCR Purification Kit (50)	50 QIAquick Spin Columns, Buffers, Collection Tubes (2 ml)	28104
QIAquick PCR Purification Kit (250)	250 QIAquick Spin Columns, Buffers, Collection Tubes (2 ml)	28106
QIAquick Gel Extraction Kit (50)	50 QIAquick Spin Columns, Buffers, Collection Tubes (2 ml)	28704
QIAquick Gel Extraction Kit (250)	250 QIAquick Spin Columns, Buffers, Collection Tubes (2 ml)	28706
QIAquick Nucleotide Removal Kit (50)	50 QIAquick Spin Columns, Buffers, Collection Tubes (2 ml)	28304
QIAquick Nucleotide Removal Kit (250)	250 QIAquick Spin Columns, Buffers, Collection Tubes (2 ml)	28306

* Routine purifications from gel slices >100 mg which contain fragments <100 bp or where the gels contain >2% agarose require additional Buffer QX1 to perform the full number of extractions.

Notes

QIAGEN International Sales



Germany QIAGEN GmbH
USA QIAGEN Inc.
Canada QIAGEN Inc.
France QIAGEN S.A.
Japan QIAGEN K.K.
Switzerland QIAGEN AG
UK QIAGEN Ltd
Australia QIAGEN Pty Ltd
ACN 072 382 944

Orders 03-892-230
Orders 800-426-8157
Orders 800-572-9613
Orders 01-60-920-920
Telephone 03-5805-7261
Orders 061-319-30-30
Orders 01293-422-9111
Orders 03-9489-3666

• Fax 02103-892-233
• Fax 800-718-2056
• Fax 800-713-5951
• Fax 01-60-920-925
• Fax 03-5805-7263
• Fax 061-319-30-33
• Fax 01293-422-922
• Fax 03-9489-3888

• Technical 02103-892-240
• Technical 800-DNA-PREP (800-362-7737)
• Technical 800-DNA-PREP (800-362-7737)
• Technical 01-60-920-930
• Technical 03-5805-7261
• Technical 061-319-30-31
• Technical 01293-422-999
• Technical 03-9489-3666

<http://www.qiagen.com>

QIAGEN International Distributors

Austria/Hungary/Slovenia

R. u. P. MARGARITELLA
Ges. m.b.H.
BIOTRADE
Breitenfurter Straße 480
1230 Wien-Rodaun
Austria
Tel: (01) 889 18 19
Fax: (01) 889 18 19 20
E-mail: BIO-TRADE@TELECOM.AT

Belgium/Luxemburg

Westburg b.v.
P.O. Box 214
3830 AE Leusden
The Netherlands
Tel: 0800-1-9815
Fax: (31) 33-4951222
E-mail: info@westburg.nl
Web site: <http://www.westburg.nl>

Brazil

Labtrade do Brazil
Av. Barão do Rego Barros, 542
V. Congonhas
Cep: 04612-041
São Paulo-Brasil
Tel: (11) 543 1455
or 0800 551321
Fax: (11) 531 3210
E-mail: labtrade@sol.com.br
Web site: <http://www.labtrade.com.br>

Central & South America

Labtrade Inc.
6157 NW 167th Street F-26
Miami, FL 33015
USA
Tel: (305) 828-3818
Fax: (305) 828-3819
E-mail: labtrade@icanet.net
Web site: <http://www.labtrade.com>

China

Gene Company Limited
Unit A, 8/F., Shell Industrial Building
12 Lee Chung Street
Chai Wan, Hong Kong, P.R.C.
Tel: (852)2896-6283
Fax: (852)2515-9371
E-mail:
Hong Kong: genehk@hkstar.com
Beijing: gene@info.yuol.cn.net
Shanghai: gene@public.sta.net.cn
Chengdu: gene@public.cd.sc.cn
Guangzhou:
gzytiao@public1.qwangzhou.gd.cn

Czech Republic

BIO-CONSULT spol. s r.o.
Bozejovická 145
142 00 Praha-Libuš
Tel: (02)4447 1239
Fax: (02)47 29 792
E-mail: bio-cons@login.cz

Denmark

KEBO Labs A/S
Roskildevej 16
2620 Albertslund
Tel: 43 86 87 88
Fax: 43 86 87 90
E-mail: cellmol@kebolab.dk
Web site: <http://www.kebolab.dk>

Egypt

Clinilab
P.O. Box 12 El-Manial
4, 160 St., El-Elehad Square
Riham Tower, El-Maadi
Cairo
Tel: 525 7212
Fax: 525 7210
E-mail: Clinilab@intouch.co

Finland

KEBO Lab Oy
Niityrinne 7
02270 Espoo
Tel: (09)804 551
Fax: (09)804 55200
E-mail: kebo@kebolab.fi

Greece

BioAnalytica S.A.
11, Laskareos Str.
11471 Athens
Tel: (01)643 61 38
Fax: (01)646 27 48
E-mail: bioanalyt@hol.gr

India

Genetix
C-88 Lower Ground Floor
Kirti Nagar
New Delhi-110 015
Tel: (011)542 1714
Fax: (011)546 7637
E-mail: genetix@nda.vsnl.net.in

Israel

Bio-Lab Ltd.
Atarot-Industrial Area
P.O.B 34038
Jerusalem 91340
Tel: 02. 584 1111
Fax: 02. 584 1110
E-mail: bio_lab@netvision.net.il
Web site: <http://www.bio-lab.co.il>

Italy

Genenca (M-Medical srl)
Via Pier Capponi, 57
50132 Firenze
Tel: (055) 500 1871
Fax: (055) 500 1875
E-mail: m.medical@agora.stm.it
Web site:
<http://www.agora.stm.it/m.medical>

Korea

IRS Laboratories, Inc.
SongBuk P.O. Box 61
Seoul, 136-600
Tel: (02) 924-86 97
Fax: (02) 924-86 96
E-mail: inslab@nowmuri.nowcom.co.kr

Malaysia

Research Biolabs Sdn Bhd
79A Jalan SS15/4C
Subang Jaya
47500 Selangor
Tel: (03)7312099
Fax: (03)7318526
E-mail: BICLABS@TM.NET.MY

Mexico

Quimica Valaner
Jalapa 77 Col. Roma
P.O. Box 24-432
06700 México, D.F.
México
Tel: 525-5725
Fax: 525-5625
E-mail:
qvanelaner@infosel.net.mx/QIAGEN

The Netherlands

Westburg b.v.
P.O. Box 214
3830 AE Leusden
Tel: (033)4950094
Fax: (033)4951222
E-mail: info@westburg.nl
Web site: <http://www.westburg.nl>

New Zealand

Biolab Scientific Ltd.
244 Bush Road
Albany, Auckland
Tel: (09)980-6700
or 0800-933-966
(09)980-6788
E-mail: info@biolab.co.nz
Web site: <http://www.biolab.co.nz>

Norway

KEBO Lab AS
Postboks 45, Kallbakken
0901 Oslo
Tel: 22 90 00 00
Fax: 22 90 00 40
E-mail: kebo@kebolab.no

Poland

Syngen Biotech
ul. Legnicka 62a
54-204 Wrocław
Tel: (071) 351 41 06
or 351 54 20
Fax: (071) 351 04 88
E-mail: syngen@infonet.wroc.pl
Web site: <http://www.syngen.com.pl>

Portugal

IZASA PORTUGAL, LDA
Rua Cordeiro Ferreira, 9
1700 Lisboa
Tel: (1)-751 6000
Fax: (1)-759 9529

Singapore

Research Biolabs Pte Ltd.
29 Lucky Crescent
Singapore 467742
Tel: 445 7927
Fax: 448 3966
E-mail:
BICLABS@SINGNET.COM.SG

Slovak Republic

BIO-CONSULT Slovakia s.r.o.
Ruzovna dolina 6
SK-824 70 Bratislava
Tel/Fax: (07) 50221 336
E-mail: bio-cons@login.cz

South Africa

Southern Cross Biotechnology (Pty) Ltd
P.O. Box 23681
Claremont 7735
Cape Town
Tel: (021) 671-5166
Fax: (021) 671-7734
E-mail: info@scb.co.za
Web site: <http://www.scb.co.za>

Spain

IZASA, S.A.
Aragón, 90
08015 Barcelona
Tel: (93) 902.20.30.90
Fax: (93) 902.22.33.66

Sweden

KEBO Lab
Fagerstogatan 18A
16394 Spånga
Tel: (08) 621 34 00
Fax: (08) 621 34 70
E-mail: cellmol@kebolab.se

Taiwan

TAIGEN Bioscience Corporation
P.F. No. 306, Section 4, Chen-Der
Road
Taipei
Taiwan, R.O.C.
Tel: (02) 2880 2913
Fax: (02) 2880 2916
E-mail: taigen@ms110.hinet.net

Thailand

Theera Trading Co. Ltd.
64 Charan Sanit Wong Road
(Charan 13) Bangkokkaiy
Bangkok 10600
Tel: (02) 412-5672
Fax: (02) 412-3244
E-mail: theeratrad@samart.co.th

QIAGEN Importers

Saudi Arabia

Abdulla Fouad Co. Ltd.
Medical Supplies Division
P.O. Box 237, Damman 31411
Tel: (03) 8324400
Fax: (03) 8346174

Turkey

DATEKS
4. Cadde 141/2
06510 Emek-Ankara
Tel: (312) 212 70 97
Fax: (312) 212 07 82

All other countries

QIAGEN GmbH, Germany

